


## Product data sheet - ReadyCell

The Sepax protocol software is an integral part of the Sepax system and provides a fully-automated method to separate the cellular product into its components, allowing a safe, reproducible and user independent procedure.

ReadyCell is a protocol designed for concentration of the Total Nucleated Cell (TNC) fraction from various cellular products into a small final volume of 7-20 ml.

The ReadyCell protocol is particularly suitable for bone marrow volume reduction, also in point of care scenarios.

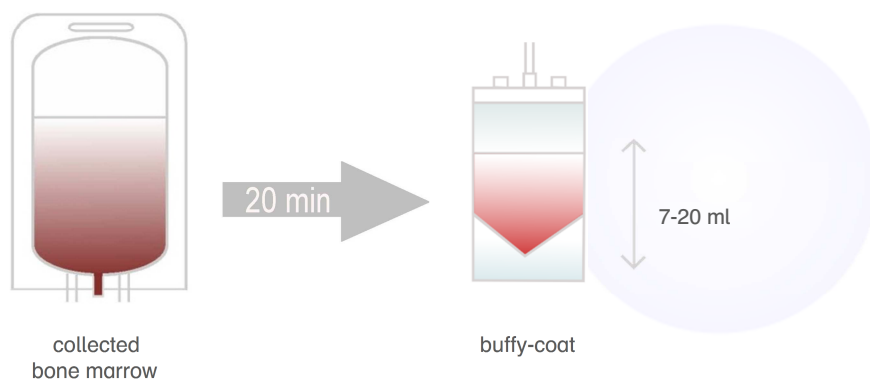

ReadyCell features	
Processed initial volume	30 to 220 ml
Selectable final volume	7 to 20 ml
Processing time	~15 minutes
TNC concentration factor	Between 3 and 10

### Specifics


#### Step-by-step work flow

1. After bone marrow aspiration and filtration, connection of product bag to single-use kit
2. Installation of kit on Sepax device
3. Automated procedure
4. End of procedure: concentrated cells are easily collected from single-use kit vial, ready for injection

#### Sepax Device settings:

- Traceability IDs: Possibility to use the traceability kit of the Sepax (barcode reader + printer) in order to trace specific identification criteria.
- Summary print-out: Possibility to print a report at the end of the procedure with the traceability kit of the Sepax (barcode reader + printer).
- Protocol available in: English, French, German, Spanish and Italian.

Associated single-use kit: CS-470.1 (refer to dedicated product datasheet)

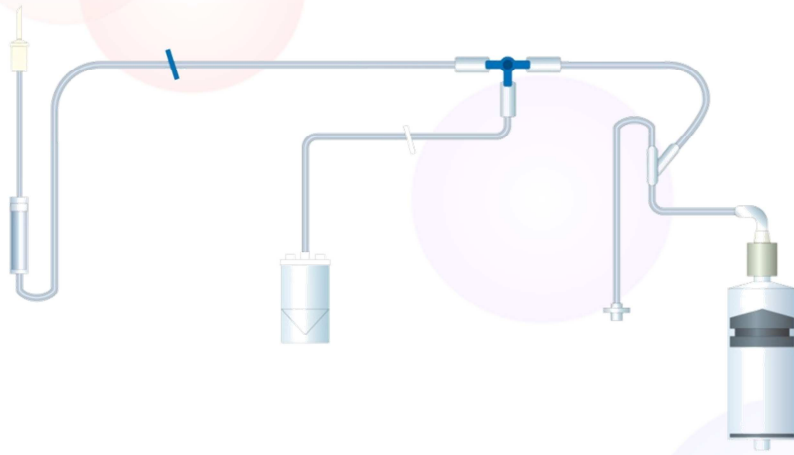

**Single-use kit x protocol software matrix**

Kit/ protocol	ReadyCell	GVR	NeatCell	CordWash	SmartWash
CS-470.1	X				
CS-490.1		X			
CS-600.1				X	X
CS-900.2			X		

**Order information:**

ReadyCell  
Packaging

Product #14302  
box including protocol software on USB key and user guide for installation

Head Office  
**Biosafe Group SA**  
Eysins, Switzerland  
T: +41 22 365 27 27  
E: info@biosafe.ch

North America Office  
**Biosafe America, Inc.**  
Houston, USA  
T: +1 713 936 2615  
E: info@biosafeamerica.com

Asia office  
**Biosafe Asia-Pacific Ltd.**  
Hong Kong  
T: +852 2956 7500  
E : info@biosafeasiapacific.com